NEW SMART® REAGENT SYSTEMS

TEST FACTOR	REAGENT SYSTEM CODE NUMBER	TEST NUMBER
Alkalinity	3670-SC	001 Alkalinity T
Hardness, Total	3671-SC	047 Hardness T
pH, Phenol Red	3672-SC	076 pH T
Cyanuric Acid	3673-SC	040 Cyanuric Acid T
Urea	3674-SC	096 Urea

The above SMART Reagent Systems can be used with the BT family of LaMotte meters and the BLE family of LaMotte meters.

The meter family identity can be found in three places:

- The splash screen when the meter is powered on.
- The firmware version number that is visible on the bottom left corner of the display when the meter is powered on. The firmware version number for a BLE meter will start with a 3. The firmware version number for a BT meter will start with a 2.
- The label on the bottom of the meter.

SMART3 BLE meter (Code 1910-01)

The calibration curves for the above reagent systems are included in the firmware for the SMART3 BLE meter. The test procedures are included in the SMART3 BLE manual. They can also be found in the SMART3 BLE manual [1910-01-MN] and as individual test procedure instructions at www.lamotte.com.

SMART3 BT meter (Code 1910)

The calibration curves for the above reagent systems are NOT included in the firmware for the SMART3 BT meter. To update the firmware/software contact Software Support at 1-800-344-3100 option 2.

The test procedures are NOT included in the SMART3 BT manual. The test procedures can be found in an updated SMART3 BT manual [1910-MN-V3] and as individual test procedure instructions at www.lamotte.com.

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800-344-3100 · +1 410-778-3100 (Outside U.S.A.) · Fax: +1 410-778-6394