

# SMART3 BLE- QUICK START GUIDE

## LAMOTTE PRE-PROGRAMMED TESTS

1.	Press and briefly hold  to turn meter on. Main Menu will appear.
2.	Press  to select <b>Testing Menu</b> .
3.	Press  to select <b>All Tests</b> menu.
4.	Press  or  to scroll to the desired test.
5.	Press  to select the test.
6.	Insert the blank. Close the lid. Press  to scan the blank.
7.	Insert the reacted sample. Close the lid. Press  to scan the sample. The result will appear on the screen.

## ADD OR DELETE A TEST IN A SEQUENCE

1.	Press and briefly hold  to turn meter on. Main Menu will appear.
2.	Press  or  to the <b>Setup Menu</b> . Press  .
3.	Press  to select <b>Edit Sequences</b> .
4.	Press  or  to scroll to the desired sequence and press  .

### Add A Test

5.	To add a test before or after an existing test, press  or  to scroll to the existing test.
6.	Press  to select the existing test. Press  or  to scroll to <b>Insert Before</b> or <b>Insert After</b> .
7.	Press  to select the option. The <b>All Tests</b> menu will appear.
8.	Press  or  to scroll to the test that will be added to the sequence.
9.	Press  to select the test. The test will be added to the sequence.

### Delete A Test

5.	To delete a test, press  or  to scroll to the test in the sequence.
6.	Press  to select the test. Press  or  to scroll to <b>Delete</b> .
7.	Press  to select <b>Delete</b> . The test will be deleted from the sequence.

# CREATE A USER TEST

1. Press and briefly hold  to turn meter on. Main Menu will appear.
2. Press  or  to the **Setup Menu**. Press .
3. Press  or  to scroll to the **Edit User Test**. Press .
4. Press  or  to scroll to the desired test. Press .

## Name The Test

5. Press  to select **Name The Test**.
6. Press ,  and  to select the characters in the test name.
7. Press  after the last character has been selected.

## Select Vial/WL

8. Press  or  to select Vial/WL. Press .
9. Press  or  to scroll to the channel with the desired wavelength and vial size combination. Press .

## Enter 2 Point Calibration

10. Press  or  to scroll to **STD Calibration**. Press .
11. Press ,  and  to select the characters in the low concentration followed by the high concentration. Press .
12. Insert the blank. Press . Repeat with reacted low and high concentration standards. Press .

## Select Decimal Places

13. Press  or  to scroll to **Decimal Places**. Press .
14. Press ,  and  to select the desired number of decimal places. Press .

## Select Units

15. Press  or  to scroll to **Select Units**. Press .
16. Press  or  to select the desired units. Press . The meter will return to the **Edit Test** menu. Press  three times to return to the **Main Menu** to begin testing.

NOTE: Use the AC wall adapter to rapidly charge the meter. Reinsert the USB port plug after charging.

WARNING: Only use the USB cable [1720-01] that is supplied with the kit. Make no substitutions.



802 Washington Ave · Chestertown · Maryland · 21620 · USA  
800-344-3100 · +1 410-778-3100 [Outside U.S.A.] · Fax: 410-778-6394 · [www.lamotte.com](http://www.lamotte.com)

© 2023 LaMotte Company · 1910-BLE-QG · 07.13.2023