




Nitrate
Nitroso
Nitrate

25 Test Strips
Tiras Reactivas
Bandelettes d'Analyse

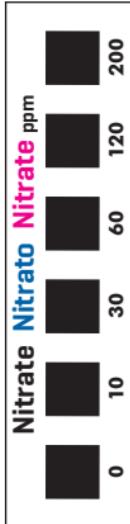

Dip pad into sample and immerse for 2 seconds. Remove strip with pad face up.

DO NOT SHAKE OFF EXCESS WATER.

In shade, hold pad level for 60 seconds and read the strip.

• Keep wet fingers out of the bottle.

• Close cap tightly after removing strip.

• Store in a cool, dry place.

Sumerja la tira en la muestra durante 2 segundos. Retire la con el lado coloreado hacia arriba. NO SACUDA EL EXCESO DE AGUA. Mantenga la tira horizontal a la sombra por 60 segundos y haga la lectura.

• No toque el envase con los dedos húmedos.

• Cierra bien el envase luego de sacar la tira.

• Guarde el envase en un lugar frío, seco y oscuro.

Immerger pour 2 secondes dans l'échantillon d'eau.

Retirer la bande face vers le haut. NE PAS SECOUER. A l'ombre, garder le niveau de la bande pour 60 secondes et lire.

• Ne pas introduire les doigts trempés dans la bouteille.

• Refermer le capuchon.

• Entreposer dans un endroit frais, sec et ombragé.

Lamotte Company
802 Washington Avenue, Chestertown, MD 21620 USA
www.lamotte.com/insta

Code 3012-G



6 37395 45337 3