

QUANTITY	CONTENTS	CODE
2 x 30 mL	*Phosphate Acid Reagent	*V-6282-G
5 g	Phosphate Reducing Reagent	V-6283-C
2	Test Tubes, 2.5-5-6.5-10 mL, glass, w/cap	0844
1	Pipet, 1.0 mL, plastic	0354
1	Spoon, 0.1 g, plastic	0699
1	Low Range Comparator Viewer	1102
1	Phosphate Low Range Comparator Bar 0.0-2.0 ppm	3122-01



*Reagent is a potential health hazard. **READ SDS:** lamotte.com
Emergency information:
Chem-Tel USA 1-800-255-3924
Int'l, call collect, 813-248-0585



To order individual reagents or test kit components, use the specified code number.


Warning! This set contains chemicals that may be harmful if misused. Read cautions on individual containers carefully.
Not to be used by children except under adult supervision.

NOTES:

- This test determines levels of orthophosphates only.
- This test should be run on clear samples only. Filter the sample if necessary.
- Best results are obtained when solution temperatures are 23-25°C.

USE OF THE LOW RANGE COMPARATOR

The Low Range Comparator allows extremely faint colors to be matched to color standards by viewing the reaction down the length of the tube increasing the path length by ten times compared to comparators that view the reaction across the diameter of the tube.

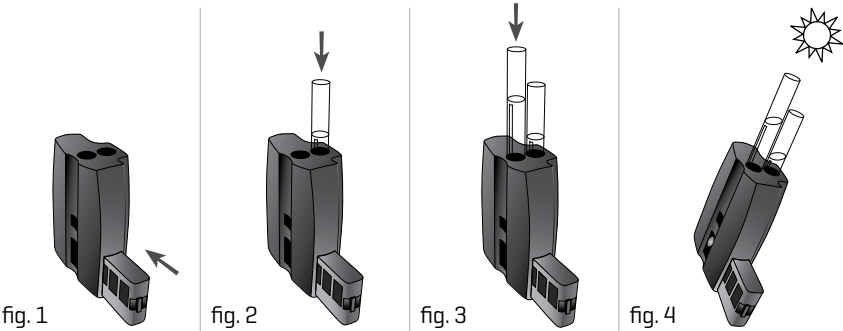

1. Slide the Low Range Comparator Bar into the Low Range Comparator Viewer [fig.1].
2. Fill a test tube to the 10 mL line with untreated sample water. Insert it in the rear hole on the top of the Low Range Comparator [fig.2].
3. Fill a test tube with sample water. Follow the test procedure. Remove the cap. Insert the tube in the front hole on the top of the Low Range Comparator [fig.3].
4. Position the comparator so that light shines down through the test tubes [fig.4]. Tilt the comparator until the color standards and sample are illuminated. Match the color of the reaction to the color standards. Read the result from the Low Range Comparator Bar.

PROCEDURE

1. Slide the Phosphate Low Range Comparator Bar [3122-01] into Low Range Comparator Viewer [1102]. Fill one test tube [0844] to the 10 mL line with sample water. Place in Low Range Comparator.
2. Fill test tube [0844] to 10 mL line with sample water.
3. Use 1.0 mL pipet [0354] to add 1.0 mL of *Phosphate Acid Reagent [V-6282]. Cap and mix.
4. Use 0.1 g spoon [0699] to add one level measure of *Phosphate Reducing Reagent [V-6283]. Cap and mix until dissolved. Wait 5 minutes.
5. Remove cap from test tube. Place tube in the Low Range Comparator. Match sample color to a color standard. Record as ppm Orthophosphate.