



## PHOSPHATE KIT

OCTA-SLIDE 2, 10-80 ppm

CODE 4401-02

QUANTITY	CONTENTS	CODE
30 mL	*VM Phosphate Reagent	*4410-G
1	Pipet, 1.0 mL, plastic	0354
2	Test Tubes, 5 mL, plastic, w/caps	0106
1	Octa-Slide 2 Viewer	1101
1	Phosphate Octa-Slide 2 Bar, 10-80 ppm	4412-01

\*WARNING: Reagents marked with an \* are considered to be potential health hazards. To view or print a Safety Data Sheet (SDS) for these reagents go to [www.lamotte.com](http://www.lamotte.com). Search the four digit reagent code number listed on the reagent label, in the contents list or in the test procedures. Omit any letter that follows or precedes the four digit code number. For example, if the code is 4450WT-H, search 4450. To obtain a printed copy, contact LaMotte by e-mail, phone or fax.


Emergency information for all LaMotte reagents is available from Chem-Tel: (US, 1-800-255-3924) (International, call collect, 813-248-0585)

To order individual reagents or test kit components, use the specified code number.

Note: The test should be run on clear samples only. Filter the samples if necessary.

Warning! This set contains chemicals that may be harmful if misused. Read cautions on individual containers carefully. Not to be used by children except under adult supervision.

### USE OF THE OCTA-SLIDE 2 VIEWER

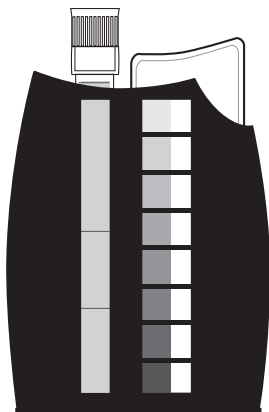

The Octa-Slide 2 Viewer should be held so non-direct light enters through the back of the Viewer. Insert the Octa-Slide 2 Bar into the Viewer. Insert the reacted sample into the top of the Viewer. Match the color of the reaction to the color standards.

## PROCEDURE

1



Insert Phosphate Octa-Slide 2 Bar (4412-01) into the Octa-Slide 2 Viewer (1101).

2




Fill a test tube (0106) to the 5 mL line with sample water.

3



Use the 1.0 mL pipet (0354) to add 1.0 mL of \*VM Phosphate Reagent (4410).

4

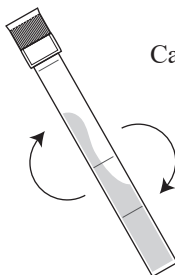

Cap and mix.

5




Wait 5 minutes.

6



Insert test tube into Octa-Slide 2 Viewer (1101).

7

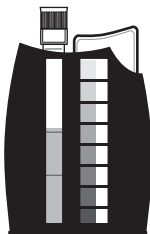

Match sample color to a color standard. Record as ppm Phosphate.

### LaMOTTE COMPANY

Helping People Solve Analytical Challenges

802 Washington Ave · Chestertown · Maryland · 21620 · USA  
800-344-3100 · 410-778-3100 [Outside U.S.A.] · Fax 410-778-6394  
Visit us on the web at [www.lamotte.com](http://www.lamotte.com)