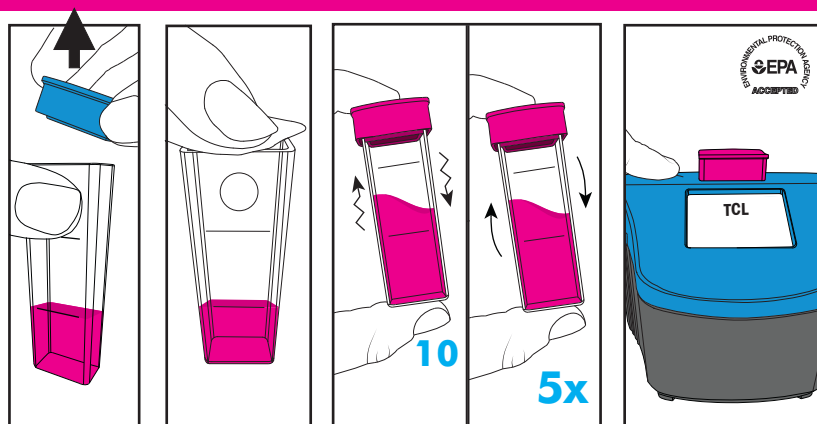
3 TCL

DPD 3 IG 6197A-J/6197A-CN-J

1. Remove cap from reacted FCL tube.
2. Add 1 **DPD 3 IG** Tablet to the same tube.
3. Cap tube with **pink** cap. Shake for 10 seconds. Invert slowly 5 times.
4. Insert tube into the ColorQ 2x.
5. Press button to read "TCL" Total Chlorine.
6. Press button to return to "Blank" or hold button until ☺ is displayed. Remove tube.

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0 - 8 ppm FCL and TCL results EPA Accepted. See www.lamotte.com for details.

INTERFERENCES: bromine, chlorine dioxide, iodine, oxidized manganese and chromium, and ozone.

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- Read the instructions before testing.
- Press button to turn meter ON and OFF ☺
- Quickly press button twice to skip a test.
- Cap the tubes securely to prevent leaks.
- Do not interchange color-coded caps.
- Dry the tubes before placing them in the chamber.
- Rinse tablet crusher after each use. Brush and rinse tubes, caps and tablet crusher after testing.
- Allow the tubes to dry before returning them to the case.
- The tube specified in the test procedure must be used.
- Use only LaMotte reagents. Cap securely.
- If unit should auto-off during testing, rinse and fill a tube. Repeat Blank and return to testing.
- ☺ = Cap tube to eliminate stray light.

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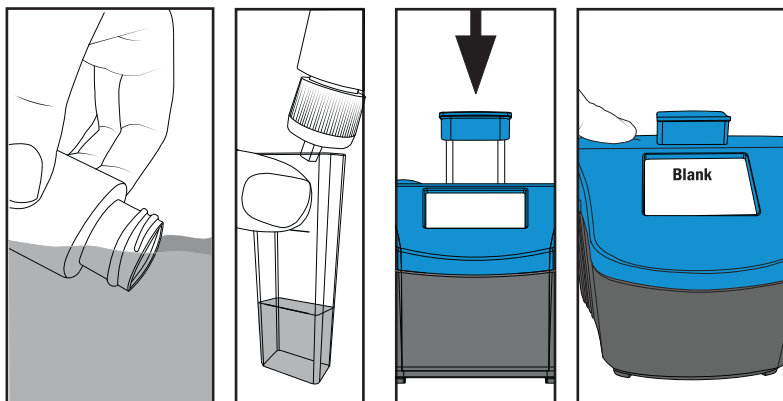
1 Blank

1. Fill sample bottle (**0689**) with water sample. Replace cap.
2. Fill clean tube (**0205**) to the 5 mL line with water sample. Cap tube with **blue** cap.
3. Insert tube into the ColorQ 2x.
4. Press button to turn meter on . When "Blank" appears press button to "Blank" the meter, and go to Free Chlorine (FCL).

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WARNING! This set contains chemicals that may be harmful if misused. Read cautions on individual containers carefully
Not to be used by children except under adult supervision.

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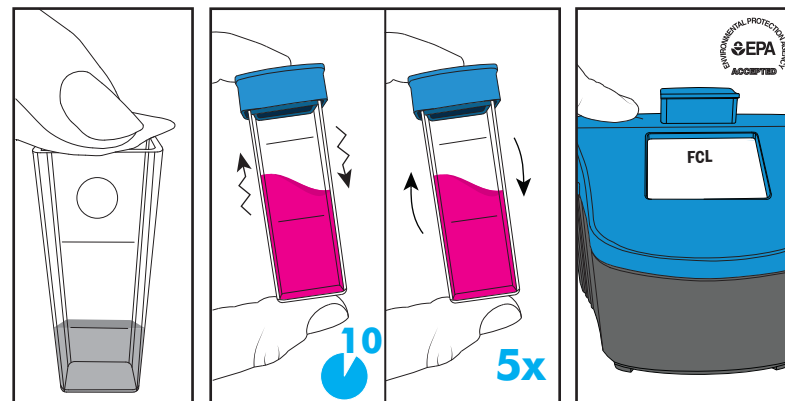
ColorQ 2x Range/Distancia/Gamme
0-10 ppm Chlorine/Chloro/Chlore**2 FCL**

DPD 1 IG 6903A-J/6903A-CN-J

1. Add 1 **DPD 1 IG** Tablet to same tube.
2. Cap tube with **blue** cap. Shake for 10 seconds. Invert slowly 5 times to mix.
3. Insert tube into the ColorQ 2x.
4. Press button to read "FCL" Free Chlorine.
5. Press button to go to "TCL" Total Chlorine. Remove tube.

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*Reagent is a potential health hazard.

READ SDS: lamotte.com **Emergency information:**
Chem-Tel USA 1-800-255-3924 Int'l, call collect, 813-248-0585

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©2023 LaMotte Company