

Code 2984LR

Peroxide Low Range  **ANALYTIC**

25 Test Strips

Insta-Test[®]

Peroxide ppm		Code 2984LR					
0	1	3	10	30	50		

Using a cup-size sample,
immerse strip for **2 seconds**.
Remove strip with **pad face up**.
DO NOT SHAKE OFF
EXCESS WATER. With pad
still facing up, immediately
compare to color chart.
Discard strip after
30 seconds.

Keep wet
fingers out of
the bottle.
Cap bottle
tightly.
Store in a cool,
dry place.

Made in
USA

LaMotte Company
802 Washington Ave., Chestertown, MD 21620 USA
410-778-3100 • www.lamotte.com

