

Code 3002

High Range
Chlorine Dioxide
50 Test Strips




Insta-Test[®] ANALYTIC

High Range Chlorine Dioxide



Using a cup-size sample,
immerse strip for **2 seconds**.
Remove strip with **pad face up. DO NOT SHAKE OFF EXCESS WATER**. With pad still facing up, **wait 10 seconds and read the strip**. Discard strip after 30 seconds.

Keep wet fingers out of the bottle.
Cap bottle tightly.
Store in a cool, dry place.
In concentrations greater than 500 ppm the pad will be a mottled color of golden-brown and black.

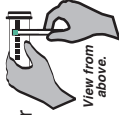

LaMotte Company
Chestertown, MD 21620
www.lamotte.com

View from
above.

