


WIDE RANGE pH TEST KIT
Code 3353-01 | Octa-Slide 2, pH 5.0-10.0






QUANTITY	CONTENTS	CODE
30 mL	*Wide Range pH Indicator	*2218-G
2	Test Tubes, 2.5-5-10 mL, plastic, w/caps	0106
1	Wide Range pH Octa-Slide 2 Bar, 5.0-10.0	3483-01
1	Octa-Slide 2 Viewer	1101

To order individual reagents or test kit components, use the specified code number.



*Reagent is a potential health hazard. **READ SDS:** lamotte.com

Emergency information:
Chem-Tel USA 1-800-255-3924
Int'l, call collect, 813-248-0585






USE OF THE
OCTA-SLIDE 2 VIEWER

Warning! This set contains chemicals that may be harmful if misused.
Read cautions on individual containers carefully.
Not to be used by children except under adult supervision.

PROCEDURE


1.

Insert Wide Range pH Octa-Slide 2 Bar [3483-01] into the Octa-Slide 2 Viewer [1101]




2.

Fill a test tube [0106] to the 10 mL line with sample water.




3.

Add 8 drops of *Wide Range pH Indicator [2218].




4.

Cap and mix.

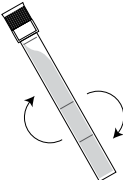

5.

Insert test tube into Octa-Slide 2 Viewer [1101]. Hold the viewer so that non-direct light enters through the back.



6.

Match sample color to a color standard. Record as pH.

