

MOLYBDENUM LOW RANGE KIT



Code 3628-01 | Test Strips

QUANTITY	CONTENTS	CODE
50	Molybdenum Test Strips	3025
30 mL	*Molybdenum Buffer	*7646-G
2	Test Tube, with caps	0106
1	Pipet, 0.5mL, plastic	0353
1	Color Chart	3628-CC

To order individual reagents or test kit components, use the specified code number.



*Reagent is a potential health hazard. **READ SDS:** [lamotte.com](https://www.lamotte.com)
Emergency information:
Chem-Tel USA 1-800-255-3924
Int'l, call collect, 813-248-0585



PROCEDURE

1.
Fill a test tube [0106]
to the 5 mL line with
sample water.

2.
Use 0.5 mL pipet [0353] to
add 0.5 mL of *Molybdenum
Buffer [7646].

3.
Remove one Molybdenum Test
Strip [3025] from vial. Gently
bend strip approximately in
half [do not fold] with pad
facing inward and visible
above the other end of
the strip. Place strip
inside test tube cap.

4.
Cap and invert gently
5 times.

5.
Remove cap.
Do not remove
strip from cap.
Wait 30 seconds.

6.
Using cap as a handle,
compare test strip to
color chart [3628-CC].
Read the result as ppm
Molybdenum [Mo].